

FROM 48 to 1, and ANYTHING in BETWEEN

Our new human panel: larger and customizable

The MILLIPLEX® Human Cytokine/Chemokine/Growth Factor Panel A combines tests for 48 individual immune factors that previously have not been together in a single panel, while also allowing you choose any combination of analytes, or select a 48-plex or a 38-plex premixed kit.

MILLIPLEX® Human Cytokine/Chemokine/Growth Factor Panel A

(Cat. Nos. HCYTA-60K*; 38-plex premixed bead kits: HCYTA-60K-PX38, HCYTA-60K-PXBK38; 48-plex premixed bead kits: HCYTA-60K-PX48, HCYTA-60K-PXBK48)


*Also available for customizing your premix!

Analytes: See table below

Samples: Serum, plasma, and cell/tissue culture supernatant

Standard Curves: Broad and fixed ranges for lot-to-lot consistency, optimized for each analyte

Standard Curves: 48 Assays

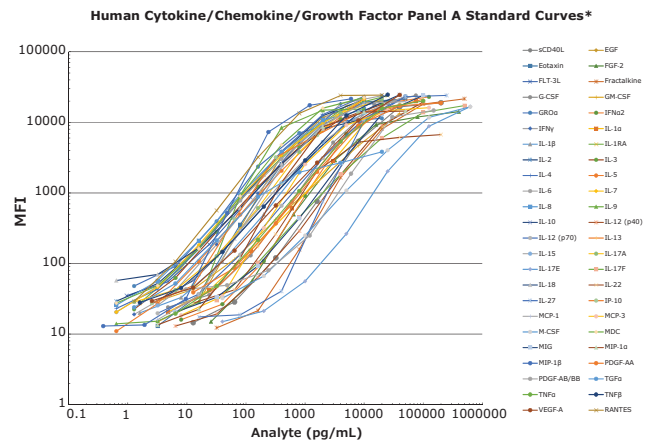

*Standard curves for all analytes were performed in MXHSM-A serum matrix except RANTES, which was in L-AB assay buffer. The kit was run using the HCYTA-60K overnight assay protocol.

Built with the trusted MILLIPLEX® assay specifications to guarantee consistent performance over time

Analyte	Standard Curve Range (pg/mL)	Sensitivity: minDC (pg/mL)*
sCD40L	13 – 200,000	5.65
EGF°	3 – 50,000	3.20
Eotaxin°	3 – 50,000	3.08
FGF-2	26 – 400,000	22.30
FLT-3L	0.96 – 15,000	0.84
Fractalkine	32 – 500,000	29.75
G-CSF°	4.8 – 75,000	3.76
GM-CSF°	2.6 – 40,000	1.55
GRO α	1.3 – 20,000	1.05
IFN α 2°	8 – 125,000	6.56
IFN γ °	1.3 – 20,000	0.86
IL-1 α °	4.8 – 75,000	2.27
IL-1 β °	1.6 – 25,000	0.52
IL-1RA°	1.6 – 25,000	1.29
IL-2°	0.64 – 10,000	0.28
IL-3°	1.3 – 20,000	0.28
IL-4°	0.64 – 10,000	0.20
IL-5°	0.64 – 10,000	0.17
IL-6°	0.64 – 10,000	0.14
IL-7°	0.64 – 10,000	0.14
IL-8°	0.64 – 10,000	0.52
IL-9	0.64 – 10,000	3.05
IL-10°	2.6 – 40,000	0.91
IL-12 (p40)°	6.4 – 100,000	3.24

Analyte	Standard Curve Range (pg/mL)	Sensitivity: minDC (pg/mL)*
IL-12 (p70)°	3 – 50,000	0.88
IL-13°	6.4 – 100,000	2.58
IL-15°	3 – 50,000	0.74
IL-17A°	1.3 – 20,000	0.71
IL-17E/IL-25°	40 – 625,000	19.77
IL-17F°	32 – 500,000	28.63
IL-18°	0.64 – 10,000	0.53
IL-22°	13 – 200,000	12.68
IL-27	16 – 250,000	50.78
IP-10/CXCL10°	2.6 – 40,000	2.13
MCP-1°	3 – 50,000	3.05
MCP-3	8 – 125,000	8.61
M-CSF°	40 – 625,000	31.95
MDC	0.64 – 10,000	0.42
MIG°	6.4 – 100,000	3.98
MIP-1 α °	3 – 50,000	3.82
MIP-1 β °	0.38 – 6,000	0.37
PDGF-AA°	13 – 200,000	10.33
PDGF-AB/BB°	9.6 – 150,000	16.39
RANTES°	1.3 – 20,000	1.58
TGF α	1.3 – 20,000	0.97
TNF α °	6.4 – 100,000	5.39
TNF β °	1.6 – 25,000	0.80
VEGF-A°	2.6 – 40,000	0.98


*Assay sensitivities (minDC, or minimum detectable concentrations) calculated using the overnight assay protocol with n = 8 assays, with MILLIPLEX® Analyst 5.1 software, view our protocol for more information.

°Contained in the 38-plex premixed kit.

so Many analytes!

Same ready-to-use kit format with a straightforward protocol


Testing Disease Samples: Sepsis vs. Normal

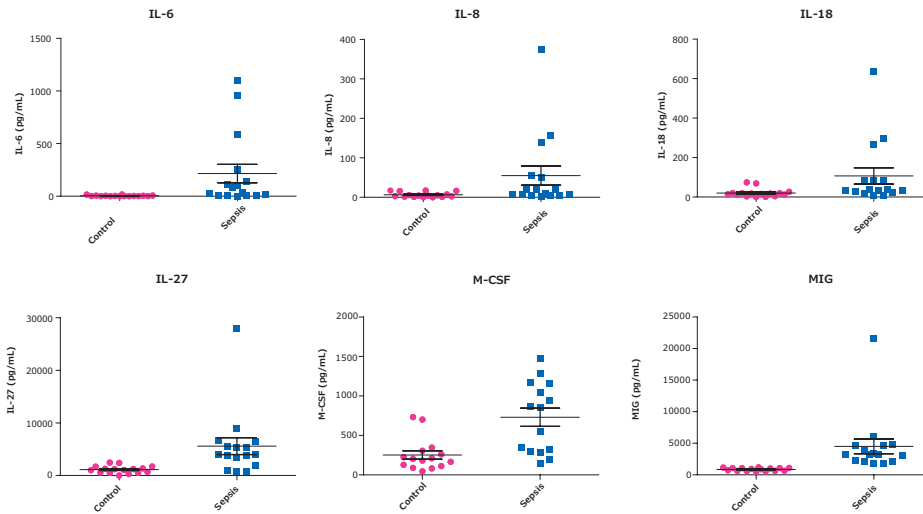

Healthy control serum/plasma samples (obtained from BioIVT) and sepsis patient serum/plasma samples (obtained from BioIVT, Discovery, and BioCemed) were tested neat (25 μ L/well) in the HCYTA-60K panel.

Healthy control serum/plasma samples, N=20. Sepsis serum/plasma samples, N=16.

Testing Stimulated PBMCs

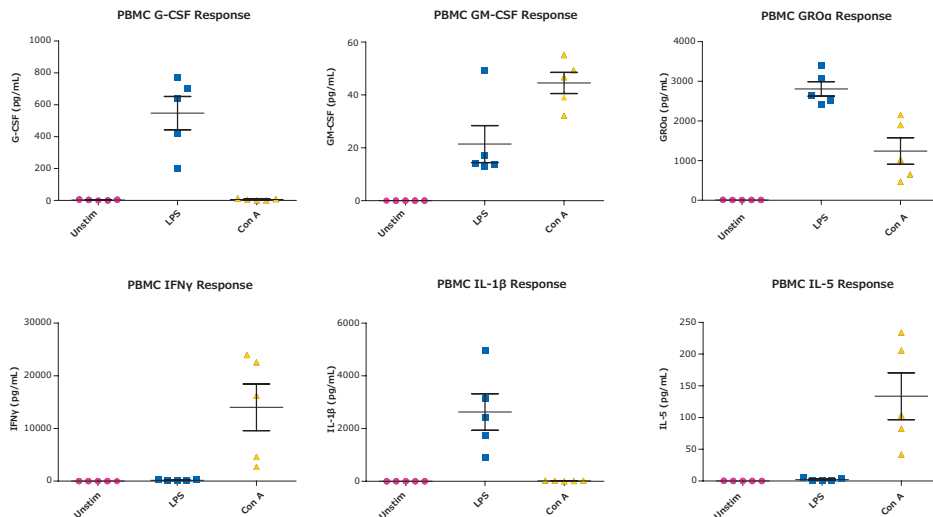

Human PBMC samples (obtained from BioIVT) in RPMI containing 10% FBS and 1% Penicillin/ Streptomycin at 10^6 cells/mL were incubated with either 1 μ g/mL LPS or Con A (Sigma-Aldrich) for 48 hours at 37 $^{\circ}$ C, after which cell-free supernatants were collected and tested with the HCYTA-60K panel.

MILLIPLEX[®] kits are manufactured in facilities which are ISO 9001-2015 compliant and are for research use only (RUO), not for use in clinical or for medical diagnostic purposes.

To view our large portfolio of MILLIPLEX[®] assays for Luminex[®] technology, visit

SigmaAldrich.com/milliplex

To place an order or receive technical assistance

Order/Customer Service: SigmaAldrich.com/order
 Technical Service: SigmaAldrich.com/techservice
 Safety-related Information: SigmaAldrich.com/safetycenter

SigmaAldrich.com

Merck KGaA
 Frankfurter Strasse 250
 64293 Darmstadt, Germany

© 2019 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. Merck, the vibrant M and Millipore are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

Lit. No. MK_FL4916EN Ver. 1.0 2019 - 25765 10/2019



Distributed by Abacus dx

1800 ABACUS (AUS) 0800 222 170 (NZ) | info@abacusdx.com | www.abacusdx.com

abacus dx