

Light the way to YOUR NEXT DISCOVERY

SMC® Neurofilament Light Chain (NF-L) Assay

Neurofilament Light Chain (NF-L) is one of the most highly studied neuroscience biomarkers in both academic and clinical research. To support your neuroscience research needs, we offer the SMC® Human NF-L High Sensitivity Immunoassay Kit (Cat. No. 03-0202-00). Single Molecule Counting (SMC®) kits using ultrasensitive immunoassay technology and the SMCxPRO® instrument platform, enable accurate, precise, and reproducible measurement of NF-L in human serum, plasma, and cerebrospinal fluid (CSF) samples.

Every SMC® kit undergoes a comprehensive development and manufacturing process to meet our in-depth verification criteria:

- Spike and recovery
- Dilutional linearity
- Parallelism in normal and diseased samples
- Intra- and inter-assay precision
- Temperature and freeze/thaw stability testing

The SMC® NF-L Assay Features Comprehensive Verification to Ensure Your Success

Standard Curve Range	0.43 pg/mL – 1,500 pg/mL
Lower Limit of Quantification (LLOQ) ¹	0.87 pg/mL
Inter-Assay %CV	9.60%
Sample Volume (µL) (Serum; K2 EDTA Plasma; CSF)	50; 50; 25
Sample Values: Mean (Range) pg/mL (Serum; K2 EDTA Plasma; CSF)	8.1 (1.9-21.9); 2.7 (1.4-3.8); 97.5 (85.4-180.3)
Spike Recovery ² : Mean % (Serum; K2 EDTA Plasma; CSF)	93; 94; 123
Dilution Linearity ³ : Mean % (Serum; K2 EDTA Plasma; CSF)	110; 106; N/A
Parallelism ⁴ : Mean % (Serum; K2 EDTA Plasma; CSF)	N/A; N/A; 99

Assay Verification Characteristics for the SMC® NF-L Assay. Thorough characterization is performed for all SMC® high sensitivity immunoassays to ensure reliable measurements in human serum, K2 EDTA plasma, and cerebrospinal fluid.

1. Lower limit of quantitation (LLOQ): lowest point on standard curve with CV <20% and accuracy within 20% of expected values

2. Spike recovery: expressed as % accuracy of measurement of recombinant standard analyte in each sample matrix serially-diluted starting from the recommended dilution factor

3. Dilution linearity: expressed as % accuracy of measurement of endogenously expressed analyte in each reported sample matrix serially-diluted starting from the recommended dilution factor

4. Parallelism: expressed as % accuracy of measurements of endogenously expressed analyte serially-diluted starting from the recommended dilution factor

Kit Ordering Information

Product Description	Cat. No.
SMC® Human NF-L High Sensitivity Immunoassay Kit	03-0202-00

For Research Use Only. Not For Use In Diagnostic Procedures.



Accurate Sample Measurements

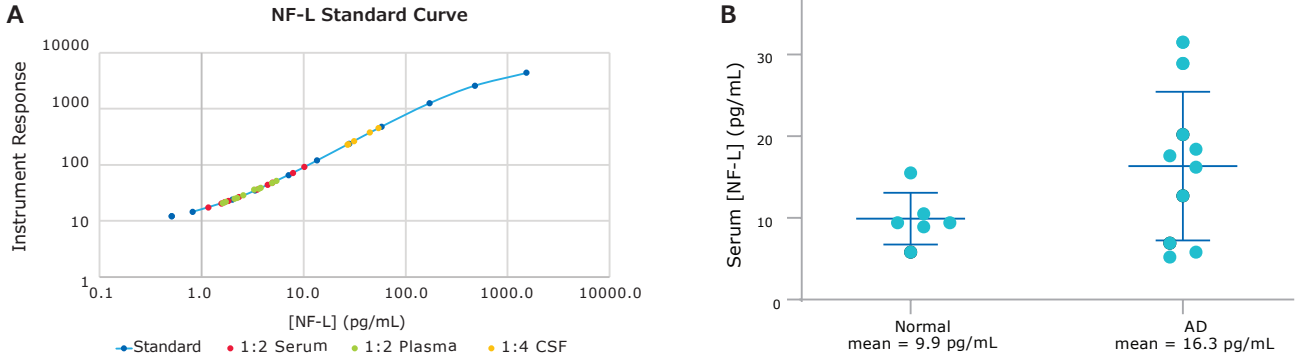


Figure 1. Sample data from the SMC® NF-L assay. Alzheimer's disease (AD) and (normal/non-AD) human samples were commercially sourced and run in the SMC® NF-L assay according to protocol. **(A)** Serum (n=10, red), K2 EDTA plasma (n=10, green), and CSF (n=5, yellow) samples accurately fall within the standard curve range (blue). **(B)** NF-L serum concentrations of AD (n=10) show nearly double that of normal (n=6) samples.

Competitor Assay Correlation

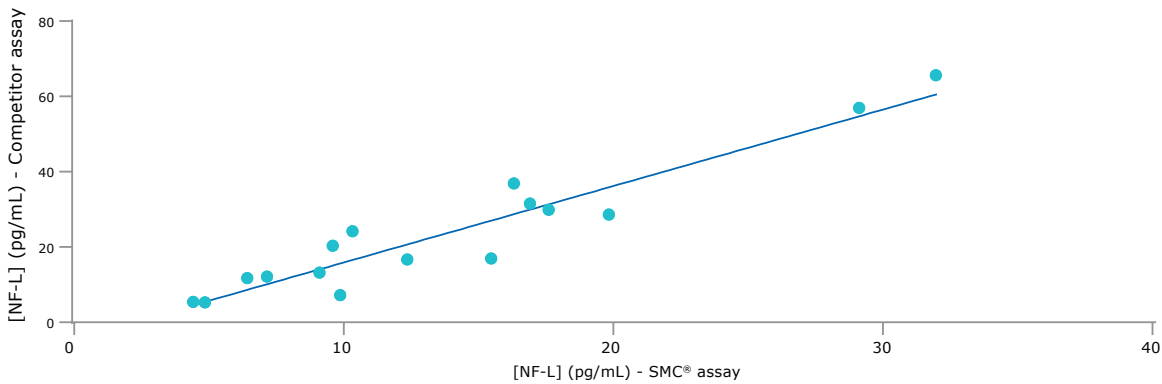


Figure 2. The SMC® NF-L assay correlates with the leading competitor kit. Human serum samples (n=6 normal/non-AD and n=10 AD) were commercially sourced and run according to protocols in the SMC® NF-L assay and in a NF-L kit offered by a leading competitor. NF-L concentration values were interpolated from each assay standard curve and plotted as shown. Good correlation ($R^2=0.92$) exists between the two assays.

Discover more at
SigmaAldrich.com/smc-neuro



To place an order or receive technical assistance
 Order/Customer Service: SigmaAldrich.com/order
 Technical Service: SigmaAldrich.com/techservice
 Safety-related Information: SigmaAldrich.com/safetycenter

Merck KGaA
 Frankfurter Strasse 250
 64293 Darmstadt, Germany

SigmaAldrich.com

We have built a unique collection of life science brands with unrivalled experience in supporting your scientific advancements.

Millipore® Sigma-Aldrich® Supelco® Milli-Q® SAFC® BioReliance®

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. Merck, the vibrant M, BioReliance, Millipore, Milli-Q, SAFC, Sigma-Aldrich, Supelco, and SMCxPRO are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.
 MK_FL9863EN Ver. 1.0 44246 11/2022



Distributed by Abacus dx

1800 ABACUS (AUS) 0800 222 170 (NZ) | info@abacusdx.com | www.abacusdx.com

abacus dx