

Cell Connect™

The Cell Connect™ is designed to streamline the addition of hematopoietic stem cells and freezing solution into up to four CryoStore™ freezing bags to prepare for cryopreservation.

It is designed to be sterile docked to both your stem cells and freezing media for closed system transfer of products, potentially eliminating the need for a clean room.




The benefits of using Cell Connect™ include:

- Ease of use saves on training time needed for laboratory personnel
- Efficiency - make 6 connections and you are done
- Closed system reduces contamination issues
- Needle-free processing reduces potential for injury
- PVC tubing allows for sterile docking

Cell Connect™

Example of use:

Close all pinch clamps. Attach cells to spike (A) and attach freeze media to spike (B). Attach up to four CryoStore™ freezing bags via male luers (E). Seal off any unused tubes to reduce cell loss. Open pinch clamps to the desired pathway, transfer cells and solutions. A rocker or oscillator is recommended for freeze media addition.

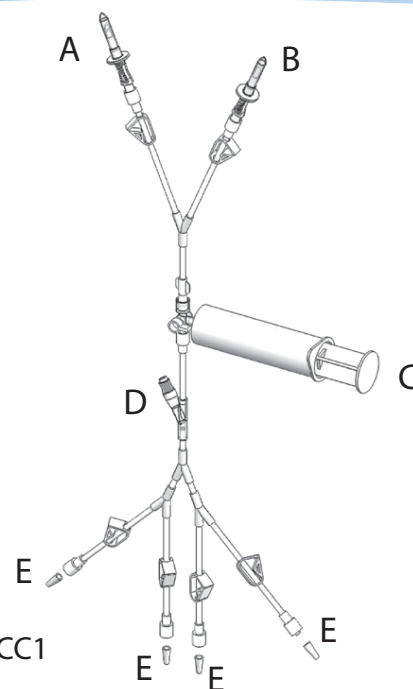

Product Code: CC1

| Product Code | Description | Stock/Non-Stock | Case Qty |
|--------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|-----------------|----------|
| CC1 | Two spikes to a 3-way stopcock, 60ml syringe and needle-free valve to four male luers with 70cm SCD tubing and pinch clamps. | Stock | 12 |
| CC2 | One spike and male luer to a 3-way stopcock, 60ml syringe and needle-free valve to four male luers with 70cm SCD tubing and pinch clamps. | Stock | 12 |
| CC35-5MT | 5 male luers with pinch clamps. Intended as an add-on to CC1 or CC2 via sterile dock to create 5 additional terminals for a total of 8 SCD connections for CryoStore™ bag filling. | Non-Stock | 12 |

Individually packed in Tyvek® pouch, supplied sterile, for single use only.



7000 Burleson Rd, Bldg D
Austin, TX USA 78744-3202
Tel: +1 (512) 474 7278
Tel: Toll Free (800) 233 9014
E-mail: sales.us@origen.com

Cell Connect™ is a registered trademark of OriGen Biomedical Inc
© 2014 OriGen Biomedical. All Rights Reserved

Distributed by Abacus dx

1800 ABACUS (AUS) 0800 222 170 (NZ) | info@abacusdx.com | www.abacusdx.com

abacus dx